

Startup and Shutdown procedure for the MA6

- **Startup [Staff Only]**
 - Turn the Tool **ON** in Coral.
 - Power Up the MA6 by turning the **POWER SWITCH ELECTRONICS** to the **ON** position.
 - Turn **ON** the Constant Intensity Controller (CIC) by pressing **ON** button on the front panel.
 - The software version is shown on the display.
 - The CIC performs a self-calibration test and displays “**ready.**”
 - Press **CP** (constant power) key. Display shows **wait**, followed by **START**.
 - Press **START** key. This will ignite the exposure lamp.
 - **LED LAMP LIFE/POWER** is flashing until lamp warming up is finished.
 - Let the lamp warm up for 15 minutes.
 - Press **CH1** to set the tool up for constant intensity.
 - Use the standard operation procedure for normal use.
- **Shutdown [Staff Only]**
 - Turn **OFF** the Constant Intensity Controller by pressing **OFF** on the front panel.
 - Power down the MA6 machine by turning the **POWER SWITCH ELECTRONICS** to **OFF** position.
 - **Shutdown** the tool on Coral.